

PeptiGel[®] Protocol:

How to Process PeptiGel[®] Cultures for Cell Staining

Before you begin, please note:

This protocol describes how to process PeptiGel[®] cultures for cell staining including histology, immunohistochemistry and immunofluorescence.

Two options, wax embedding and freezing for cryo-sectioning, are outlined here.


It's recommended to use Greiner Bio-One ThinCerts[™] for this protocol. These inserts are made with polystyrene and will dissolve during the xylene wash step.

This protocol has been written as a guide. Therefore, further assay optimization may need to be carried out for your cultures.


Protocol 1: Wax Embedding

Fixation:

1. Prewarm 10% neutral buffered formalin (NBF) solution to room temperature. Immerse the PeptiGel[®] (+/- cells) into the solution and incubate for 30 minutes at room temperature (~25°C).

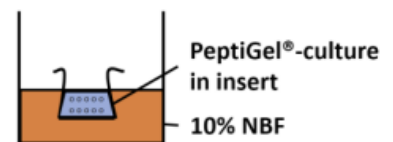

2. Replace the NBF solution with phosphate buffered saline (PBS) solution, enough to immerse the PeptiGel[®]. Store at 4°C until processing in wax. The maximum recommended storage time is 2 weeks.

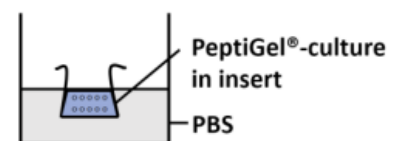

Processing in Wax:


3. Place the fixed insert containing the PeptiGel[®] cultures (+/- cells) in a cassette. Then place the cassette in a processor basket and assemble into the processor.




- Run the wax processing according to the guideline in the table below.

Reagent	Time (seconds)
50% IMS	90
70% IMS	60
99%IMS	60
99%IMS	60
99%IMS	60
99%IMS	60
99%IMS	60
Xylene	90
Xylene	90
Xylene	90
Molten-wax	90
Molten-wax	120

- Embed the sample according to your instrument's instructions.

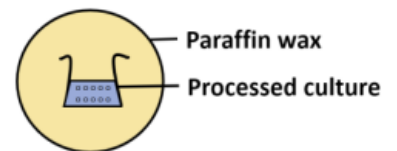



- Cryo-section according to your instrument's instructions. Individual sections of 5-7 µm thickness are recommended.

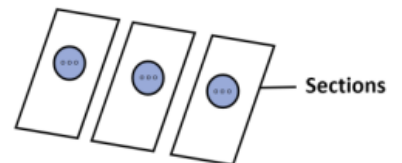


Protocol 2: Freeze for Cryo-Sectioning

Fast Freeze Using Isopentane:


- Decant liquid nitrogen (LN) into a wide mouthed vacuum flask. Place 25-50 mL of isopentane in a small thick-walled glass beaker and immerse in LN to cool. Lift out when isopentane starts to freeze around the rim of the beaker.
- Apply a drop of optimal cutting temperature (OCT) compound on to a 20 mm cork disc. Place a PeptiGel® (+/- cells) sample onto the disk and apply more OCT until the sample is covered.

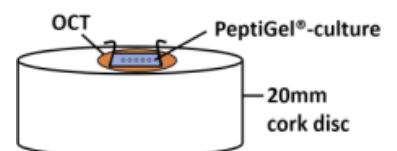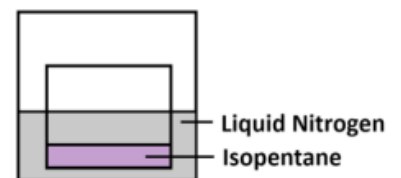

- Place PeptiGel® (+/- cells) with the disc into the isopentane for 5-15 seconds to freeze, then remove.

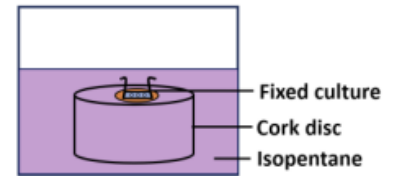


- Wrap the frozen PeptiGel® in aluminium foil and store at -80°C.

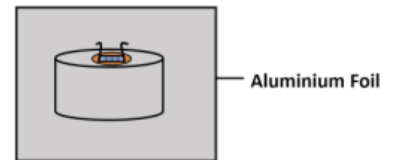

Sectioning Using Cryostat:

- Cryosection according to the instrument's instructions. Cutting sections of 7-10 µm or cutting a 20 µm thickness if culture samples lack structure, is recommended.

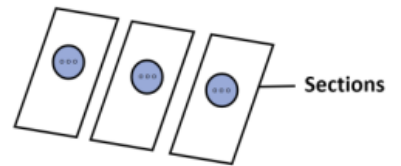

For further support, please contact our technical support team at tech@cellgs.com.



General info@cellgs.com
Technical Enquiries tech@cellgs.com
Orders order@cellgs.com

www.cellgs.com

EUROPE
Cell Guidance Systems Ltd
Maia Building
Babraham Bioscience Campus
Cambridge
CB22 3AT
United Kingdom
T +44 (0) 1223 967316
F +44 (0) 1223 750186

USA
Cell Guidance Systems LLC
Helix Center
1100 Corporate Square Drive
St. Louis
MO 63132
USA
T 760 450 4304
F 314 485 5424